

Package contents

Verify the shipping list (included with shipment) corresponds with the crate contents. Contact apricotsupport@sptlabtech.com immediately if any items are missing.

WARNING!



**Contents are heavy.
2 persons required.**

Unpacking the crate

- ① With a Phillips screwdriver, remove the 28 packaging screws from all four sides of the crate.
- ② Lift off the top of the crate and remove any boxes and accessories.
- ③ Lift the sides of the crate up and off the base and remove the tool kit.
- ④ With the 14mm open-ended wrench, remove all four bolts that secure the instrument to the bottom of the crate.
- ⑤ Remove the transit blocks by pushing in, then pull down and out.

